


Deliver medication confidently with **MediGel**[®] Sucralose

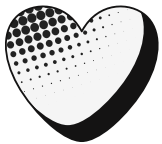

Flexible Dosing and Easy Intake

MediGel[®] Sucralose is a flavored, low-calorie water gel that masks bitter medications and promotes voluntary intake.



Improve Outcomes and Reduce Stress

The soft gel texture allows for thermoreversible mixing of drugs directly into the cup, no water bottles, gavage, or injections required.



Simplify Preparation and Administration

Simply mix your medication into the liquid gel, allow it to solidify, and serve to your animals. See the full mixing instructions on the next page.



Thermoreversible Gel Mixing Instructions




Step 1

Place the gel cup in a water bath at the liquefaction temperature and time, i.e. 60° C for 15 min.




Step 2

Remove from the water bath and dry the cup. Prepare the cup for medication by injection.

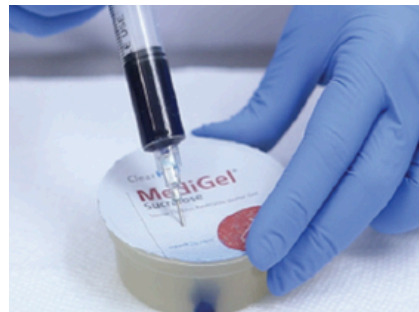


Step 3

Using a syringe with your dosage determined for a 2 oz cup of gel, carefully inject the medication.

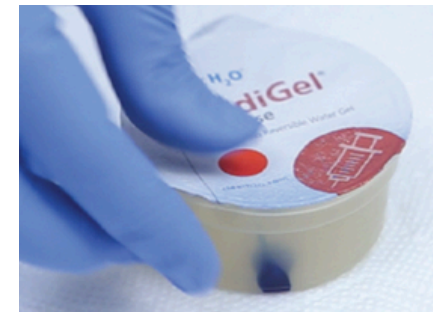

Step 4

Place a sticker over the injection site and ensure good adhesion.

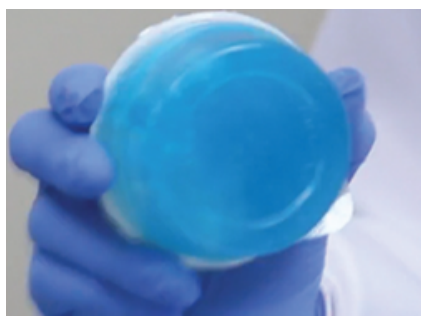

Step 5

Hand shake for a count of 15 seconds.



Step 6

Vortex for an additional count of 15 seconds.



Step 7

Date the cup. Allow the cup to cool at room temperature for at least 15 minutes.



Step 8

To accelerate solidification, place the cup in a refrigerator.